

Provitro AG

Charitéplatz 1

tel +49.30.450 578 358

sales@provitro.de www.provitro.de

Charité Campus Mitte

10117 Berlin

fax +49.30.450 578 919 www

Skeletal muscle cell growth medium, basal

Cat.-Nr.: 200 0602

contains of:

Basal media		Supplements
200 0602	500 ml Skeletal muscle cell growth medium,	-
	basal	

Maintenance of skeletal muscle cell growth medium:

Place the bottle of basal medium in the dark at 4°C to 8°C immediately after delivery.

Characteristics:

The Provitro skeletal muscle cell growth medium is a sterile liquid culture medium for culturing human skeletal muscle cells (HSKMC). The medium is delivered as a basal medium and is suitable for culturing HSKMC **after adding the optional available essential supplement kit components**. The final formulation is optimized for initial seeding of 4,000 cells / cm² up to confluence (approx. 90 %). Feeder-layer, matrix substrates or other substances are not necessary.

Stability and storage:

The supplemented skeletal muscle cell growth medium can be stored in the dark at 4°C to 8°C for up to 1 month. Do not heat the medium over 37°C or use uncontrollable sources of heat (e.g. microwave appliances). If only a part of the medium is to be used, remove this amount from the bottle and heat it.

Special note:

Do not freeze the medium. This can lead to high salt concentrations by freezing out pure water which will cause irreversible damage.

Quality control:

Provitro's skeletal muscle cell growth medium is thoroughly tested after each production. All components are tested in a stringent biological assay. Each batch is checked for HSKMC proliferating characteristics. The cells cultured in skeletal muscle cell growth medium are checked regarding their morphology, the adherence rate, the colony forming efficiency and the population doubling time.

Product specification:

The pH is set at 7.6 and osmolality at $285 \pm 10 \text{ mOsm} / \text{kg}$.

In vitro laboratory use only.

Not intended for any human or animal diagnostic or therapeutic use.