

Provitro AG Charité Campus Mitte Charitéplatz 1 10117 Berlin tel +49.30.450 578 358

fax +49.30.450 578 919

sales@provitro.de www.provitro.de

Endothelial cell proliferation medium, FCS-kit

Cat.-Nr.: 211 0001

contains of:

Basal media		Supplements	
200 0001	500 ml Endothelial cell proliferation medium,	222 1000	L-Glutamine
	basal	231 3500	FCS (foetal calf serum)
		226 0500	Heparin
		244 0500	human rec. EGF (epidermal growth factor)
		245 0250	human rec. bFGF (basic fibroblast growth factor)
		241 0025	human rec. VEGF
		242 0500	human rec. Long R3 IGF-1
		223 0005	Ascorbic acid
		224 0010	Hydrocortisone
		236 0350	Antibiotics (optional)

Maintenance of endothelial cell proliferation medium:

Place the bottle of basal medium in the dark at 4°C to 8°C immediately after delivery. Store the supplements at -20°C.

Characteristics:

The Provitro endothelial cell proliferation medium is a sterile liquid culture medium for culturing endothelial cells. The medium is delivered as a basal medium and is suitable for culturing Provitro human endothelial cells after adding the supplement kit components. The formulation is optimized for initial seeding of 4,000 cells / cm² up to confluence (approx. 90 %). Feeder-layer, matrix substrates or other substances are not necessary. **Due to the possibility of reduced proliferative activity we recommend to use the antibiotic supplement for freshly isolated cells only.**

Stability and storage:

The supplemented endothelial cell proliferation medium can be stored in the dark at 4°C to 8°C for up to 1 month. Do not heat the medium over 37°C or use uncontrollable sources of heat (e.g. microwave appliances). If only a part of the medium is to be used, remove this amount from the bottle and heat it.

Special note:

Do not freeze the medium. This can lead to high salt concentrations by freezing out pure water which will cause irreversible damage.

Quality control:

Provitro's endothelial cell proliferation medium is thoroughly tested after each production. All components are tested in a stringent biological assay. Each batch is checked for human endothelial cells proliferating characteristics. The cells cultured in endothelial cell growth medium are checked regarding their morphology, the adherence rate, the colony forming efficiency and the population doubling time.

Product specification:

The pH is set at 7.6 and osmolality at 285 \pm 10 mOsm / kg.

In vitro laboratory use only.

Not intended for any human or animal diagnostic or therapeutic use.