

Provitro AG

Charitéplatz 1

tel +49.30.450 578 358

www.provitro.de

sales@provitro.de

Charité Campus Mitte 10117 Berlin

fax +49.30.450 578 919

Supplement kit for endothelial cell proliferation medium, FCS-kit

Cat.-Nr.: 215 0001

contains of:

Basal media	Supplements
-	222 1000 L-Glutamine
	231 3500 FCS (foetal calf serum)
	226 0500 Heparin
	244 0500 human rec. EGF (epidermal growth factor)
	245 0250 human rec. bFGF (basic fibroblast growth fact
	241 0025 human rec. VEGF
	242 0500 human rec. Long R3 IGF-1
	223 0005 Ascorbic acid
	224 0010 Hydrocortisone
	236 0350 Antibiotics (optional)

Maintenance of supplement kit for endothelial cell growth medium:

The supplements are delivered on dry ice. Store the supplements at -20°C.

Characteristics:

The Provitro supplement kit for endothelial cell proliferation medium is suitable for culturing Provitro human endothelial cells after adding the supplement kit components to 500 ml basal medium. The formulation is optimized for initial seeding of 4,000 cells / cm² up to confluence (approx. 90 %). Feeder-layer, matrix substrates or other substances are not necessary. Due to the possibility of reduced proliferative activity we recommend to use the antibiotic supplement for freshly isolated cells only.

Stability and storage:

Store the **supplements** at **-20°C**. Do not use the supplements after expiring date.

Special note:

Avoid any thaw/freeze cycle.

Quality control:

Provitro's supplement kit for endothelial cell priliferation medium is thoroughly tested after each production. All components are tested in a stringent biological assay. Each batch is checked for human endothelial cells proliferating characteristics. The cells cultured in endothelial cell growth medium are checked regarding their morphology, the adherence rate, the colony forming efficiency and the population doubling time.

In vitro laboratory use only.

Not intended for any human or animal diagnostic or therapeutic use.